

File commands

Save a file with a new name	Ctrl+A
Start a new file	Ctrl+N
Open an existing data file	Ctrl+O
Print the current file	Ctrl+P
Save the current file	Ctrl+S

Movement commands

Use these keys with Meander Scan, Linear Scan, or a stereological probe:	
Next Scan Site	F2
Previous Scan Site	Ctrl+F2
These keys work in both live and acquired image modes:	
Move Left	← (left arrow key)
Move Right	→ (right arrow key)
Move Up	↑ (up arrow key)
Move Down	↓ (down arrow key)

Image commands

Acquire an image stack	Ctrl+H
Display first image in the stack	Home key
Display last image in the stack	End key
Display next image in the stack	Page Up
Display the previous image in the stack	Page Down
Delete the image stack plane	Alt+U
Display Min projection	Ctrl+Shift+M
Display Max projection	Ctrl+M
Image Montage Tools	Ctrl+I
Display all color channels	F9
Display red channel	F10
Display green channel	F11
Display blue channel	F12

Serial section commands

Next serial section	Alt+Down arrow key
Previous serial section	Alt+Up arrow key
Align serial sections for 3D reconstruction	Ctrl+F

Cursor commands

These keys control the cross-hair cursor. Use the mouse wheel with mouse wheel focus turned off to adjust the circular cursor's size.

Increase cross-hair size	Ctrl+Up arrow
Decrease cross-hair size	Ctrl+Down arrow
Make cross-hair thinner	Ctrl+Left arrow
Make cross-hair thicker	Ctrl+Right arrow

Editing commands

Copy the selection	Ctrl+C or Ctrl+Insert
Paste the contents of the Windows Clipboard	Ctrl+V or Shift+Insert
Delete the selection	Del
Undo the last edit action	Ctrl+Z

Neuron tracing commands

Hide/show the tracing	Ctrl+T
Add a node	Ctrl+Alt+A
Add a tree ending	Ctrl+Alt+E
Add a spine to a tree	Ctrl+Alt+S
Add a trifurcating node	Ctrl+Alt+T
Place a varicosity	Ctrl+Alt+V

Tracing

Select the previous contour type	Alt+B
Select the next contour type	Alt+N
Previous marker type (in <u>marker placement</u> mode)	Alt+B
Next marker type (in <u>marker placement</u> mode)	Alt+N
Close current contour	C
End open contour	Esc
Hide/show tracing	Ctrl+T

Miscellaneous

Set the Z meter to 0	Alt+S
Zoom in	Ctrl + "+"
Zoom out	Ctrl + "-"
Refresh display	F5
Toggle live video	Ctrl+L
Go to Quick Launch	Ctrl+Q

Adding a custom shortcut key

1. Right-click in any ribbon bar and select **Set custom shortcuts**.
2. In the **Customize** window, select a ribbon or **All Commands** from the **Category** drop-down list.
3. Under **Commands**, select the desired command.
4. Enter the desired shortcut key in the **Press a new shortcut key** box and click **Assign**.